


What's in the Carry Case?

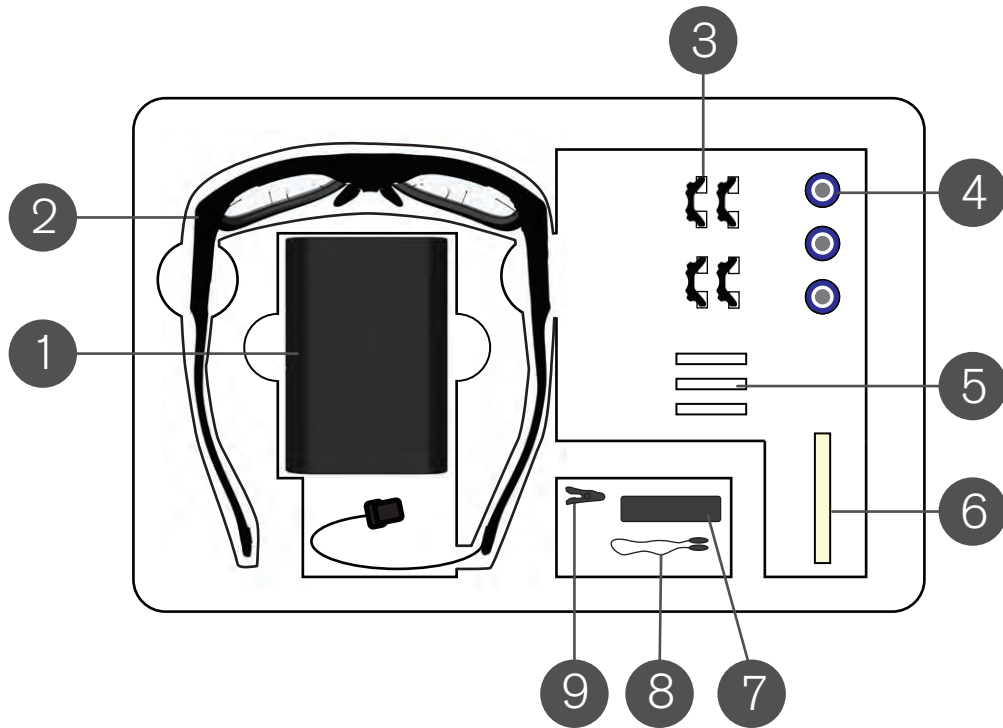

- 1 Recording unit
- 2 Head unit
- 3 4 x Nose pads
- 4 3 x Li-ion batteries
- 5 SD memory cards and sleeves
- 6 Calibration card
- 7 Microfiber cloth
- 8 Head strap
- 9 Cable clip


What's in the Accessories Box?



- 1 1 x Battery charger for Li-ion batteries
- 2 1 x USB Type-C Ethernet adapter
- 3 5 x Power cables for battery charger (EU, US, UK, AUS)
- 4 1 x USB charger for Recording unit
- 5 4 x plug adapters for USB charger (EU, US, UK, AUS)
- 6 1 x Ethernet cable (3 meters)
- 7 SD card to USB Type-A adapter
- 8 2 x Calibration cards
- 9 10 x Calibration stickers

Preparing and Connecting Tobii Pro Glasses 3

1 Prepare the recording unit

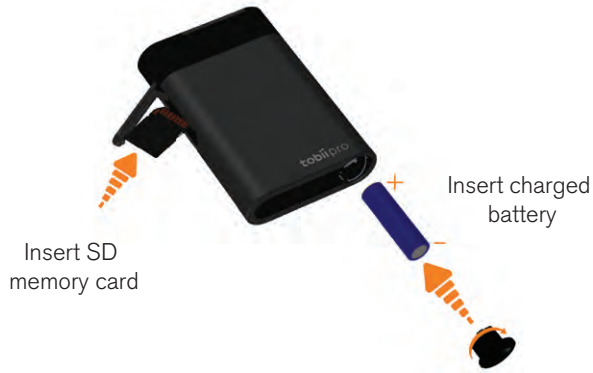

2 Install software

Follow the links below for Windows® or Android™ installation



tobiipro.com/learn-and-support/downloads-pro



play.google.com — Search for **Glasses 3**

3 Attach the nose pad (Optional)



4 Attach the head strap



5 Connect the head unit



Connect head unit cable to recording unit

6 Turn on the recording unit



Press and hold the power button for a few seconds until the power LED stops flashing

7 Connect to your device

Wireless

Find and connect to your Pro Glasses 3 via WLAN. The network name corresponds to the serial number on the back of the recording unit.

Step 1



The WLAN password for you Pro Glasses 3 is TobiiGlasses

Step 2



Wired

If using a wired connection, securely connect the Ethernet cable to the recording unit and to your computer.



Recording unit LEDs

Power Indicator LED

- (LED solid green) The recording unit is powered on. ~100-30% of battery power remaining.
- (LED solid red) The battery is running out of power. ~29-10% of battery power remaining.
- (LED blinking red) The battery is out of power. Stop any ongoing recordings immediately and replace the battery.
- (LED blinking green) The battery inside the recording unit is charging via the USB port.

SD Card Activity LED


- (LED solid green) recording unit is ready to start recording. The SD card can be removed.
- (LED solid red) There is less than enough space for 15 min. of recording.
- (LED blinking red) The SD card is running out of space, only 5 min. of recording time remains. Stop any ongoing recording immediately and replace the card.
- (LED blinking yellow) No SD card inserted, or there is something wrong with the SD card and it cannot be written to.

Connection Status LED

- (LED solid green) The recording unit is connected to the head unit
- (LED blinking yellow) The recording unit is not connected to head unit. Check the head unit cable.

Calibrate and Record with Pro Glasses 3

1 Set up recording

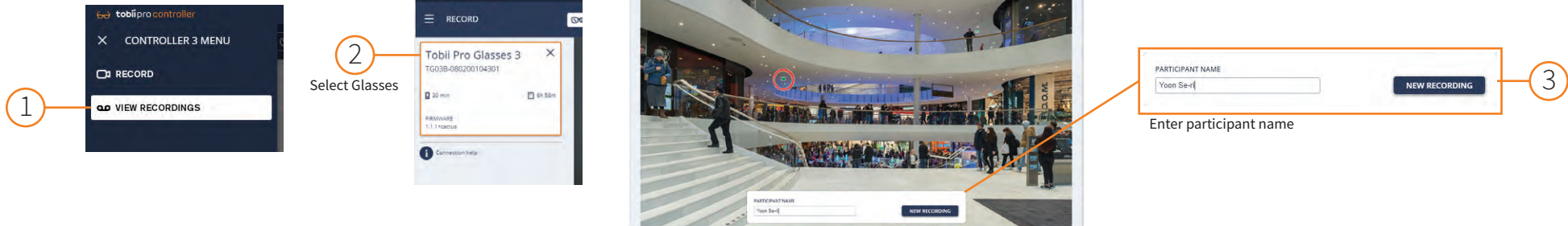


1. In the **CONTROLLER 3 MENU**, select **VIEW RECORDINGS**.

2. Select **Tobii Pro Glasses 3** (TG03B-080200104301).

3. Enter participant name in the **PARTICIPANT NAME** field and click **NEW RECORDING**.

2 Perform the calibration

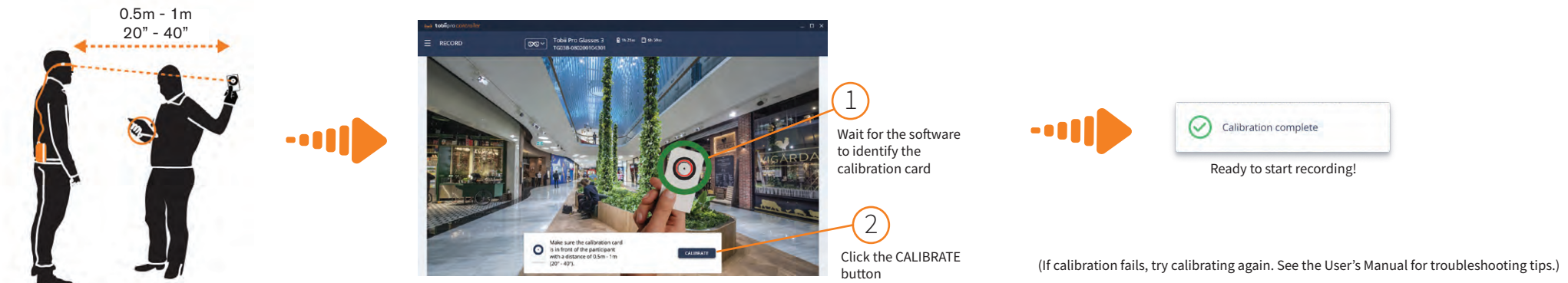

0.5m - 1m
20" - 40"


1. Wait for the software to identify the calibration card.

2. Click the **CALIBRATE** button.

Calibration complete
Ready to start recording!

(If calibration fails, try calibrating again. See the User's Manual for troubleshooting tips.)

3 Start a recording

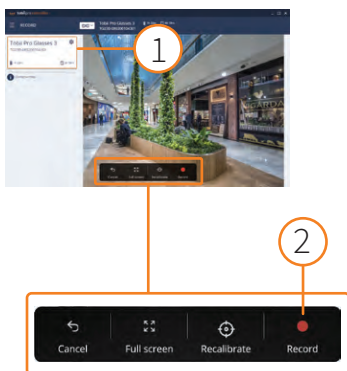

1. Start recording.

2. Record

4 Stop a recording




1. Stop Recording?

John Smith 17m 21s

2. STOP

5 Replay a recording

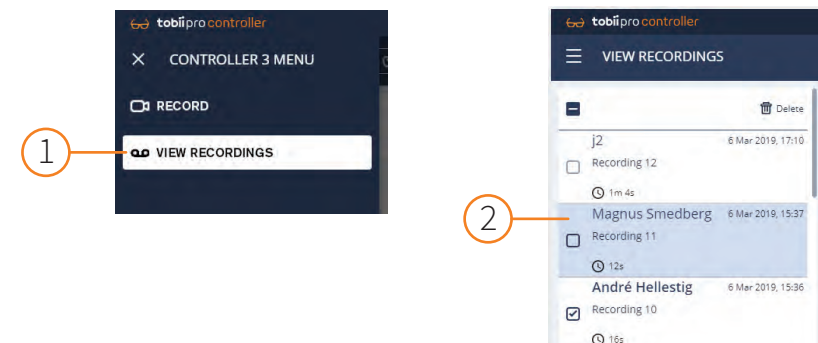

1. View recordings.

2. Select recording to replay.

Recording	Participant	Time
Recording 12	j2	6 Mar 2019, 17:10
Recording 11	Magnus Smedberg	6 Mar 2019, 15:37
Recording 10	André Hellestig	6 Mar 2019, 15:36